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**• INSTRUCTIONS ON HOW TO USE THE BUNSEN BURNER**

Appliance tested by IMQ primacontrol - test report n°55-3-PO-03

**SPECIFICATIONS :**

BURNER Mod. BUNSEN	Codice: 140 2440 01 can be used with either methane or GPL
BURNER Mod. BUNSEKUR MULTIGAS	Codice: 140 2050 10 complete with thermocouple, can be used with either methane or GPL
BURNER Mod. BUNSEKUR	Codice: 140 2050 06 complete with thermocouple, can be used with methane
BURNER Mod. BUNSEKUR	Codice: 1402050 08 complete with thermocouple, can be used with GPL

SPECIFICATIONS	G 30/G31 (GPL)	G 20 (Methane)
(Ø jets mm) only 140 2440 01	0.35	0.60
Nominal power	2.35 kW	0.70 kW
Hourly consumption	170 g/h	67 l/h
Supply pressure	1.0 bar	20 mbar

**THE BUNSEKUR MULTIGAS model is designed to work with GPL; if it is to be used with methane:**  
 Use a pair of pliers to loosen the lock nut (10)  
 Use the Allen key supplied to loosen the lock pin (8)  
 Turn the gas adjustment ring (6) and the air adjustment ring (5) anti-clockwise until the flame is blue  
 Tighten the screw (8) using the Allen key supplied.

**CONNECTING THE BUNSEN BURNER**

Only qualified technicians should connect the BUNSEN burner to the gas supply.

- Connect the rubber hose to the hose nipple on the regulator or the pipe work and tighten the hose clamp
- Connect the free end of the hose to the hose nipple on the valve of the appliance and tighten the hose clamp, make sure that the hose is not twisted.
- Keep the BUNSEN burner away from inflammable materials
- Never replace the gas bottle near sources of heat

**CONDITIONS OF USE AND CLEANING**

**LIGHTING:**

- Open the valve on the gas supply
- Open the gas valve on the BUNSEN burner by turning the knob anticlockwise, move the flame towards the burner, we recommend that a long match is used to prevent burning yourself (for BUNSEKUR models, press the safety valve prior to bringing the match near the burner and keep pressed for 20/30 seconds before releasing it).
- Use the valve on the appliance to adjust the flame according to your needs.
- After use close the valve on the gas bottle and then the valve on the burner.
- When in use, ensure that the air-intake is not blocked with rags or other objects

- If there are any gas leaks, close the valve on the gas bottle immediately
- Always use protective gloves when handling particularly hot objects
- Never move the BUNSEN burner when in use

**MAINTENANCE:**

- Check that the air-intake tube is free of dirt; Use a brush to clean if dirty.
- Periodically replace the flexible hose if it has passed the expiry date printed on the hose itself
- Periodically clean the holes in the burner using a stiff brush
- Do not move the appliance when in use
- Close the valve on the gas bottle after use.

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
Pos.	Description	Material
1	Thermocouple	Copper
2	Nozzle	Chrome plated brass
3	Thermocouple support	304 Stainless steel
4	Column	Nickel plated brass
5	Air adjustment ring	Nickel plated brass
6	Gas adjustment ring	Nickel plated brass
7	Gas adjustment needle	Brass
8	Grain	Stainless steel
9	Relief valve	Brass
10	Locknut	Nickel plated brass
11	Valve	Nickel plated brass
12	Hose nipple	Brass
13	Base	Chrome plated steel
14	Nut CH 13	Chrome plated steel

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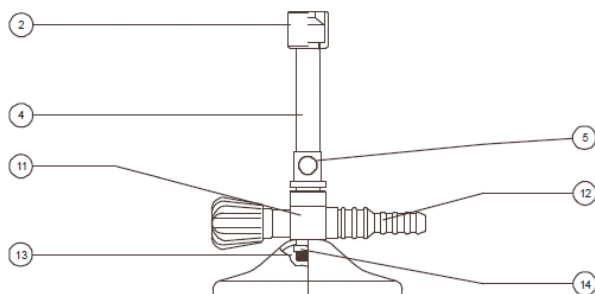Pos.	Description	Material
2	Nozzle	Chrome plated brass
4	Column	Nickel plated brass
5	Air adjustment ring	Nickel plated brass
11	Valve	Nickel plated brass
12	Hose nipple	Brass
13	Base	Chrome plated steel
14	Nut CH 13	Chrome plated steel

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