



LiquiSampler | sample collector for liquids



LiquiSampler for liquids, according to DIN EN ISO 15528. Made of ultra pure and chemically inert PTFE/FEP or PP for contamination-free sampling. Sampling from open and closed containers, barrels, tanks, silos, water courses. Suitable for all commercially available barrels and containers with opening of at least 32 mm (1.26 in.) in diameter.

Easy cleaning - all surfaces are free of pores and crevices, preventing accumulation of dirt. Only round screw threads proven in foodstuffs hygiene are used. Disassembly and cleaning are extremely simple.

- ▶ Practical one-handed use at the push of a button
- ▶ With scraping ring for marking of depth and cleaning
- ▶ Transparent
- ▶ Available in PTFE/FEP or PP transparent
- ▶ Point sampling, all-layer sampling and bottom sampling possible

Use

All-layer sample

1. Slowly insert the LiquiSampler into the liquid with the button pressed.
2. Release the push-button, the valve closes.
3. Withdraw the lance and take out the sample.

Point sampling/bottom sample

1. Insert the LiquiSampler to the targeted point/to the ground with closed valve.
2. Press the button - the liquid flows in.
3. Release the push-button, the valve closes.
4. Pull out the LiquiSampler, take out the sample.



LiquiSampler

Material	Volume ml (oz.)	Immersion depth cm (in.)	Dia. mm (in.)	Item no.
PP, transparent	150 (5.07)	60 (23.62)	32 (1.26)	5330-1060 N4
PP, transparent	250 (8.45)	100 (39.37)	32 (1.26)	5330-1100 N4
PP, transparent	500 (16.91)	200 (78.74)	32 (1.26)	5330-1200 N4
PTFE/FEP, transparent	150 (5.07)	60 (23.62)	32 (1.26)	5330-2060 N4
PTFE/FEP, transparent	250 (8.45)	100 (39.37)	32 (1.26)	5330-2100 N4
PTFE/FEP, transparent	500 (16.91)	200 (78.74)	32 (1.26)	5330-2200 N4

Accessories

Description	Item no.
Cleaning brushes 25 mm (0.98 in.) dia., length 100 cm (39.37 in.)	5304-0100 N4